# BEAD GENIE™

## Instruction Manual

EN

ES

FR

DE

IT

**Change History**

| # | Date       | Version | Descriptions         |
|---|------------|---------|----------------------|
| 1 | 2026.02.24 | A       | New release          |
| 2 | 2026.07.30 | B       | Specification Update |
|   |            |         |                      |

## Table of Contents

|                               |   |
|-------------------------------|---|
| Introduction                  | 2 |
| Package Contents              | 2 |
| Safety Information            | 2 |
| Intended Use                  | 3 |
| Installation                  | 4 |
| Specifications                | 5 |
| Control Panel                 | 5 |
| Applications                  | 6 |
| Operation Instruction         | 6 |
| Maintenance                   | 7 |
| Technical Support Information | 7 |
| Technical Data                | 7 |
| Compliance                    | 8 |

## Introduction

This manual contains installation, operation, and maintenance instructions for OHAUS BEAD GENIE™ homogenizer. Please read the manual completely before using.

## Package Contents

- Bead Beating Homogenizer
- 4 tube holders for 12x 1.5/2.0 mL microtubes, 12x 1.5/2.0 mL microtubes (horizontal), 6x 15/5 mL tubes, and 3x 50/20 mL tubes
- Power supply
- Power cord

We recommend you retain the original packaging for 90 days in case you need to return the product for any reason to your distributor or Ohaus.

## Safety Information

Safety notes are marked with signal words and warning symbols. These show safety issues and warnings. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results.

**WARNING** For a hazardous situation with medium risk, possibly resulting in severe injuries or death if not avoided.

**CAUTION** For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or minor or medium injuries if not avoided.

**ATTENTION** For important information about the product. May lead to equipment damage if not avoided.

**NOTE** For useful information about the product.

## Warning Symbols

General hazard

Electric shock hazard

## Safety Precautions

**CAUTION:** Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

Always operate unit on a level surface for best performance and maximum safety.

**DO NOT** lift unit by its attachment.

Disconnect unit from the power supply prior to maintenance and servicing.

Spills should be removed promptly.

**DO NOT** immerse the unit for cleaning.

**DO NOT** operate the unit if it shows signs of electrical or mechanical damage.

**WARNING!** The protection provided by the unit may be impaired if used with accessories not provided or recommended by the manufacturer or used in a manner not specified by the manufacturer.

**WARNING! DO NOT** use the unit in a hazardous atmosphere or with hazardous materials for which the unit was not designed.

**CAUTION:** To avoid electrical shock, completely cut off power to the unit by disconnecting the power cord from the unit or unplugging from the wall outlet.

Earth Ground – Protective Conductor Terminal

Alternating Current

## Intended Use

This instrument is intended for use in laboratories, pharmacies, schools, businesses and light industry. It must only be used for processing materials as described in these operating instructions. Any other type of use and operation beyond the limits of technical specifications, without written consent from OHAUS, is considered as not intended. This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use. If the instrument is not used according to these operating instructions, the intended protection provided by the instrument may be impaired.

## Installation

Upon receiving the unit, check to ensure that no damage has occurred during shipment. It is important that any damage that occurred during transport is detected at the time of unpacking. If you do find such damage the carrier must be notified immediately.

After unpacking, place the unit on a level bench or table, away from explosive vapors. Ensure that the surface on which the unit is placed will withstand the typical heat produced by the unit. Do not position the equipment such that it is difficult to disconnect the power cord during use. Always place the unit on a clean and sturdy work surface.

The unit is supplied with a power cord that should be plugged into a properly grounded outlet. If the supplied power cord does not meet your needs, please use an approved power cord that suits local codes and electric supply. Replacement of the cord must be made by a qualified electrician.

### Prepare for an operation:

1. Put the tube holder onto the shaft.

2. Place tubes on holder, cap end inward

3. Put the lid on and tighten the thumbscrew.


4. Check if the tubes are firmly clamped.

Specifications

| Product                | BEAD GENIE                                                                           |
|------------------------|--------------------------------------------------------------------------------------|
| Control                | Digital                                                                              |
| Accessories Included   | Tube Holder for 12x 1.5/2.0 mL, 12x 1.5/2.0 mL (horizontal), 6x 15/5 mL, 3x 50/20 mL |
| Capacity               | 12x 1.5/2.0 mL microtubes, 6x 15/5 mL tubes, 3x 50/20 mL tubes                       |
| Motion                 | Reciprocating, 19 mm                                                                 |
| Speed Range            | 500 - 4800 strokes per minute                                                        |
| Timer                  | 1 second - 99 minutes                                                                |
| Pulse - ON/OFF Time    | 1 second - 99 minutes                                                                |
| Power                  | 15V DC, 4.65A                                                                        |
| Dimensions (L x W x H) | 165 x 122 x 330 mm / 6.5 x 4.8 x 13 inch                                             |
| Net Weight             | 6 kg / 13.3 lb.                                                                      |

Control Panel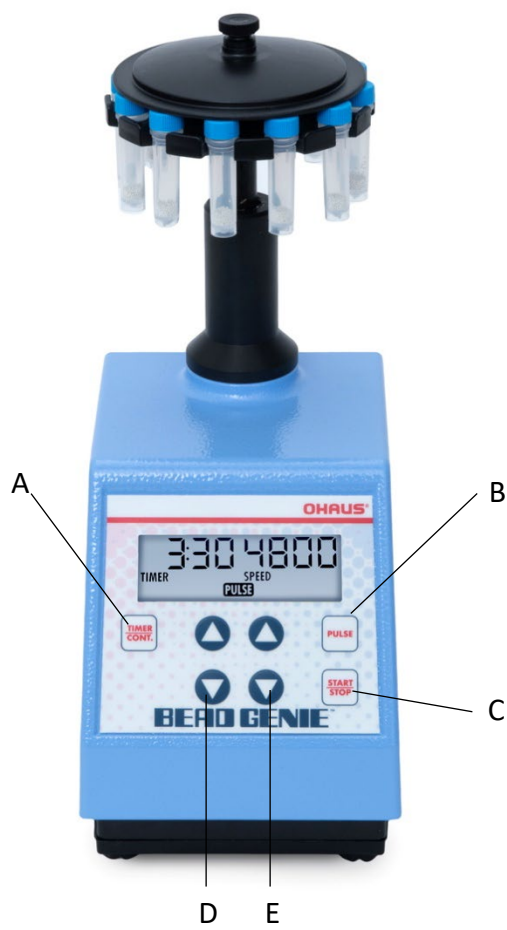

|          |                         |
|----------|-------------------------|
| <b>A</b> | Timer Button            |
| <b>B</b> | Pulse Button            |
| <b>C</b> | Start / Stop Button     |
| <b>D</b> | Timer Up / Down Buttons |
| <b>E</b> | Speed Up / Down Buttons |

## Applications

The OHAUS BEAD GENIE is a vigorous bead beating homogenizer. Applications include cell lysis, tissue homogenization, DNA/RNA extraction, protein extraction, microbial disruption, etc.

During homogenization, friction inside the sample tubes will generate heat. The BEAD GENIE is capable of being operated in a cold room or an environmental chamber. In most short runs, 3 minutes or less, it is feasible that if you refrigerate the sample tubes, beads and sample before the homogenization run, you will be able to perform the homogenization at room temperature.

The small size of the microtubes may allow sufficient heat to dissipate so that sample integrity is not impaired. The operator must make the decision based on the type of beads, the sample content and size and the length of time of the run. In any event, it is advisable to keep the tubes, beads and sample chilled before each run.

## Operation Instructions

- 1.0 - Place the BEAD GENIE on a sturdy, level work surface.
- 1.1 - Connect the power cord to a properly grounded three-pronged receptacle.
- 1.2 – Install the tube holder and tubes well by following the chapter of INSTALLATION.
- 1.3 - The BEAD GENIE is now ready for use.

## MODES OF OPERATION

2.0 - The BEAD GENIE has the following modes of operation:

### 2.1 – “TIMED”

Press the **TIMER/CONT.** button to display **TIMER** on the LCD. Press the UP/DOWN arrow button beneath the timer segment to set the desired cycle time from 1 second to 99 minutes. Next, set the speed control by pressing the UP/ DOWN arrow beneath the speed segment to any value between 500 and 4800. The setting of 500 is the least vigorous and 4800 the most vigorous. Press the **START/STOP** button to begin the bead beating action. The homogenizer will run until the set time has elapsed. To stop the cycle before the set time has elapsed, press the **START/STOP** button. To initiate a new timed cycle, press the **START/STOP** button again.

### 2.2 – “CONTINUOUS”

The “Continuous” mode allows for unattended operation that will continue indefinitely until you stop the cycle by pressing the **START/STOP** button. To initiate the “Continuous” mode, press the **TIMER/CONT.** button to display **CONTINUOUS** on the LCD. Upon pressing the **START/STOP** button, the homogenizer will start and continue to run, unattended, until you once again press the **START/STOP** button. To restart the cycle, press the **START/STOP** button again.

### 2.3 – “PULSED”

The PULSED mode will work with TIMED or CONTINUOUS modes. To initiate the PULSED mode, press the **PULSE** button to enter setting. Press the UP/DOWN arrow buttons on the left

side to **SET PULSE ON** time and press the UP/DOWN arrow buttons on the right side to **SET PULSE OFF** time. Press the **PULSE** button again to set. The word **PULSE** will be illuminated on the LCD when PULSED mode is ON.

#### 2.4 - "SLEEP"

BEAD GENIE will automatically go into a "SLEEP" mode after 10 minutes of inactivity. Press any button to "WAKE UP".

#### 2.5 - "ERROR"

When a fault is detected, BEAD GENIE will stop working and **Err** will be displayed on the LCD. Unplug to clear error and re-plug in to restart.

## Maintenance

The OHAUS BEAD GENIE homogenizer is built for long, trouble-free, dependable service. No lubrication or other technical user maintenance is required.

The unit should be given the care normally required for any electrical appliance. Avoid wetting or unnecessary exposure to fumes. Spills should be removed promptly. DO NOT use a cleaning agent or solvent on the front panel which is abrasive or harmful to plastics, nor one which is flammable. Always ensure the power is disconnected from the unit prior to any cleaning. If the unit ever requires service, contact your Ohaus representative.

## Technical Support Information

For service assistance or technical support, please visit our website at [www.ohaus.com](http://www.ohaus.com) to locate and contact the Ohaus office nearest you.

Serial Number: \_\_\_\_\_

Date of Purchase: \_\_\_\_\_

Supplier: \_\_\_\_\_

## Technical Data

### Equipment Rating

- Indoor use only
- Altitude: 0 to 2000 m
- Environmental: 0°C to 38°C (32°F to 100°F), max 95% relative humidity, non-condensing
- Overvoltage category (Installation category): II

## Compliance

Compliance to the following standards is indicated by the corresponding mark on the product.

| Mark                                                                              | Standard                                                                                                                                                                                                                                                                                                                        |
|-----------------------------------------------------------------------------------|---------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
|  | This product complies with the applicable harmonized standards of EU Directives 2011/65/EU (RoHS), (EU)2015/863, 2014/30/EU (EMC) and 2014/35/EU (LVD). The EU Declaration of Conformity is available online at <a href="http://www.ohaus.com/ce">www.ohaus.com/ce</a> .                                                        |
|  | This product complies with the EU Directive 2012/19/EU (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment.<br><br>For disposal instructions in Europe, refer to <a href="http://www.ohaus.com/weee">www.ohaus.com/weee</a> . |
|  | EN 61326-1                                                                                                                                                                                                                                                                                                                      |

### Global Notice

Warning: This is a Class A product. In a domestic environment this product may cause radio interference in which case the user may be required to take adequate measures.

### ISED Canada Compliance Statement:

CAN ICES-003(A) / NMB-003(A) FCC Notice

### FCC Compliance Statement:

This device complies with part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) This device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation.

### ISO 9001 Registration

The management system governing the production of this product is ISO 9001 certified.

**FCC Supplier Declaration of Conformity**

Unintentional Radiator per 47CFR Part B

Trade Name: OHAUS CORPORATION

Model: BEAD GENIE e-BH3...

**Party issuing Supplier's Declaration of Conformity**

Troemner LLC

201 Wolf Drive

Thorofare, NJ, 08086

USA

Phone: +1 856 686 1600

**Responsible Party – U.S. Contact Information**

Ohaus Corporation

8 Campus Drive, Ste. 105

Parsippany, NJ 07054

United States

Phone: +1 973 377 9000

Web: [www.ohaus.com](http://www.ohaus.com)

## LIMITED WARRANTY

OHAUS products are warranted against defects in materials and workmanship from the date of delivery through the duration of the warranty period. During the warranty period OHAUS will repair, or, at its option, replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to OHAUS.

This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than OHAUS. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by OHAUS Corporation. OHAUS Corporation shall not be liable for any consequential damages.

As warranty legislation differs from state to state and country to country, please contact OHAUS or your local OHAUS dealer for further details.

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P/N: 31018508 B © 2026 Ohaus Corporation, all rights reserved